

INTRODUCING

Fisherbrand™ Traceable™ 7600 Smart Wi-Fi Connected Data Logger

New features based on feedback from real-world labs make TraceableLIVE™ more effective, more efficient, and easier to set up and use. Now you can spend more time on science and less time trying to manage your monitoring system.



Traceable 7600

Smart Wi-Fi Data Logging Refrigerator/Freezer

Reimagined for Today's Clinics and Labs

This new smart Wi-Fi connected data logger offers many of the best-loved features of the previous TraceableLIVE® connected data loggers, plus much more.

- Stay connected to your critical environments no matter where you are—get email, text, and push notifications to your internet-connected device
- NEW interactive and intuitive touch screen, enhanced security, and smart probes make calibration easier
- NEW acknowledge alarms remotely via the TraceableLIVE® app
- NEW calibrated smart sensor probes with easy-grip connectors allow you to quickly swap out probes for calibration and disruption-free monitoring
- NEW open-door sensors prevent accidental temperature excursions
- NEW onboard security measures offer optional PIN protection for your settings, to prevent tampering
- NEW operates using both 2.4 and 5 GHz WiFi
- Simple, one-time setup takes only a few minutes

Choose the Traceable™ 7600 Kit That Is the Best for Your Lab

The Traceable™ 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer is available in several formats so you can pick the one that fits your lab setup. The data logger and smart sensors are also available additionally. You can add to a kit or create a system for your individual needs.



Ordering Information

Product	Cat. No.
7600 Smart Wi-Fi Data Logger	
7600 Smart Wi-Fi Data Logger	01257641
7600 Smart Wi-Fi Data Logger Kits	
7600 Smart Wi-Fi Data Logger Kit with Bottle Smart Probe and Door Sensor	01257636
7600 Smart Wi-Fi Data Logger Kit with Bullet Smart Probe and Door Sensor	01257635
7600 Smart Wi-Fi Data Logger Kit with Vaccine Bottle Smart Probe and Door Sensor	01257634
Calibrated Smart Probes and Door Sensor	
Calibrated Temperature Bottle Smart Probe	01257640
Calibrated Temperature Bullet Smart Probe	01257639
Calibrated Temperature Vaccine Bottle Smart Probe	01257638
Door Sensor	01257637

Calibration Just Got Easier with Smart Sensors



Keep the Data Logger and Swap the Sensor for More Efficient Recalibration

The Traceable™ 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer comes equipped with smart sensor probes and open-door sensors. They're called smart because each sensor is calibrated instead of the data logger.

When it's time to recalibrate, the data logger stays put, and you can swap out the sensor that's expired for one that is ready to go. It's quick, it's easy, and it poses little to no disruption to your lab's sample monitoring activity.




3 Easy Steps to Recalibrate

Recalibrating is as easy as 1, 2, 3. The visual guide below will help you have your Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer back up and running in no time.

STEP 1

Check if calibration is near expiration.

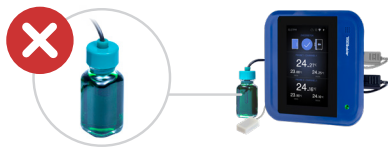

STEP 2

Purchase new probes.



STEP 3

Switch probes and you are ready to go.



Note: If you maintain a second set of probes, they can be returned to InnoCal for calibration.

Traceable® smart sensor probes and open-door sensors only work with the Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer and do not operate with other data loggers. Also, noncalibrated probes or sensors are not compatible with the Traceable® 7600 Smart Wi-Fi Data Logging Refrigerator/Freezer Thermometer.

TraceableLIVE™ Monitors Your Samples 24/7 So You Don't Have To






Know What's Happening in Your Refrigerator or Freezer, No Matter Where You Are

TraceableLIVE™ is a cloud-based subscription software solution for connected data logging devices that sends alerts to your Wi-Fi connected device when you need to take action.

- Alerts, emails, and push notifications sent to your smartphone or Apple Watch,* and voice call escalations* to your phone, for peace of mind that your samples are safe
- Wi-Fi cloud-connected, so your data is available whenever you need it
- Regular updates are automatically pushed to your system, keeping it maintenance free, up to date, and secure
- Quick to set up and easy to use, keeping you focused on what matters most
- Add multiple users and multiple devices for robust lab monitoring
- Fully compliant—meets 21 CFR Part 11 North American compliance and Annex 11 for Europe and ISO 17025 accredited calibration
- The TraceableLIVE® app is kept up to date with current technological and security developments by a team of skilled programmers and professionals

Traceable® products have an individually numbered Traceable Certificate that assures accuracy from our A2LA-accredited ISO/IEC 17025:2017 (1750.01) calibration laboratory. It indicates traceability of measurements to the SI units through NIST or other recognized national measurement institutes (NMI) that are signatories to the CIPM Mutual Recognition Agreement. Each product goes through an intense vetting process before it can carry the Traceable name, saving you both time and money by not having to calibrate separately.

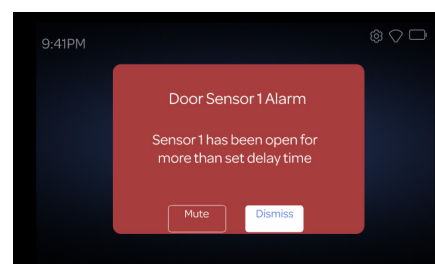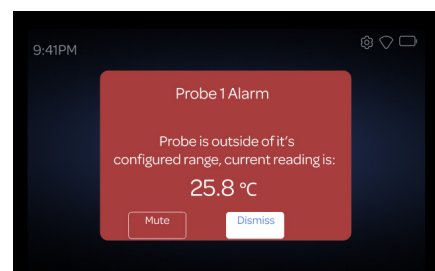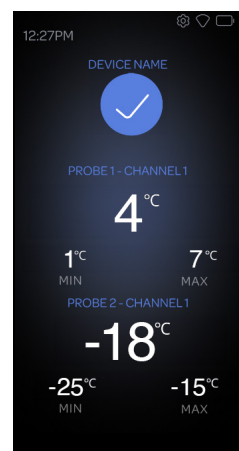

TraceableLIVE™ Subscription Plans

Purchase a 1 year or 2 year subscription plan with an activation code. TraceableLIVE® service comes in three levels. See details in the table below:

Account Information

Free Account	Standard Account	Premium Account
Try TraceableLIVE for free. Get Premium benefits for 30 days.	Equip your team to protect your sensitive items and meet most national and local regulations.**	“Future-proof” your vaccine and sample monitoring for ever-changing compliance requirements.
E-mail and push alarms	E-mail and push alarms	E-mail and push alarms
1 user only	Unlimited users	Unlimited users
1 month data storage	Unlimited data storage	Unlimited data storage
Up to 3 devices can be added to the account	No limit to devices that can be added to account (each device must have subscription)	No limit to devices that can be added to account (each device must have subscription)
Calibration due and other notifications	Calibration due and other notifications	Apple Watch compatibility
–	Pre-loaded templates set up for monitoring Spex® chemical products	SMS text notifications included
–	–	Voice notification escalation
–	–	Ideal for high importance escalations
–	–	Create your own monitoring and notification templates
–	–	Unlimited templates
–	–	Pre-loaded templates set up for monitoring Spex chemical products
–	–	21 CFR Part 11 compliance**

**Meets 21 CFR Part 11 North American Compliance and Annex 11 for Europe.

Ordering Information

Product	Cat. No.
1 Year Standard Subscription Code	01257635
2 Year Standard Subscription Code	01257634
1 Year Premium Subscription Code	01257639
2 Year Premium Subscription Code	01257638



Find out more at fishersci.com

Distributed by Fisher Scientific. Contact us today:

In the United States

Order online: fishersci.com

Call customer service: 1-800-766-7000

In Canada

Order online: fishersci.ca

Call customer service: 1-800-234-7437

