



microman® E


NEW Positive-Displacement Pipette
for viscous, volatile, and problem liquids

Full Range of Volumes

Model	Part Number	Volume Range	Price
M10E	FD10001G	1 - 10µL	\$360
M25E	FD10002G	3 - 25µL	\$360
M50E	FD10003G	20 - 50µL	\$360
M100E	FD10004G	10 - 100µL	\$360
M250E	FD10005G	50 - 250µL	\$360
M1000E	FD10006G	100 - 1000µL	\$360

Full Line of Sterilized Capillary Piston Tips

Model	Part Number	Tips per Box	Tips per Box	Price
CP10ST	F148413G	1 - 10µL	192	\$91
CP10ST	F148313G	1 - 10µL	960	\$389
CP25ST	F148712G	3 - 25µL	576	\$263
CP50ST	F148713G	20 - 50µL	576	\$263
CP100ST	F148415G	10 - 100µL	192	\$91
CP100ST	F148315G	10 - 100µL	960	\$389
CP250ST	F148714G	50 - 250µL	576	\$263
CP1000ST	F148180G	100 - 1000µL	182	\$91

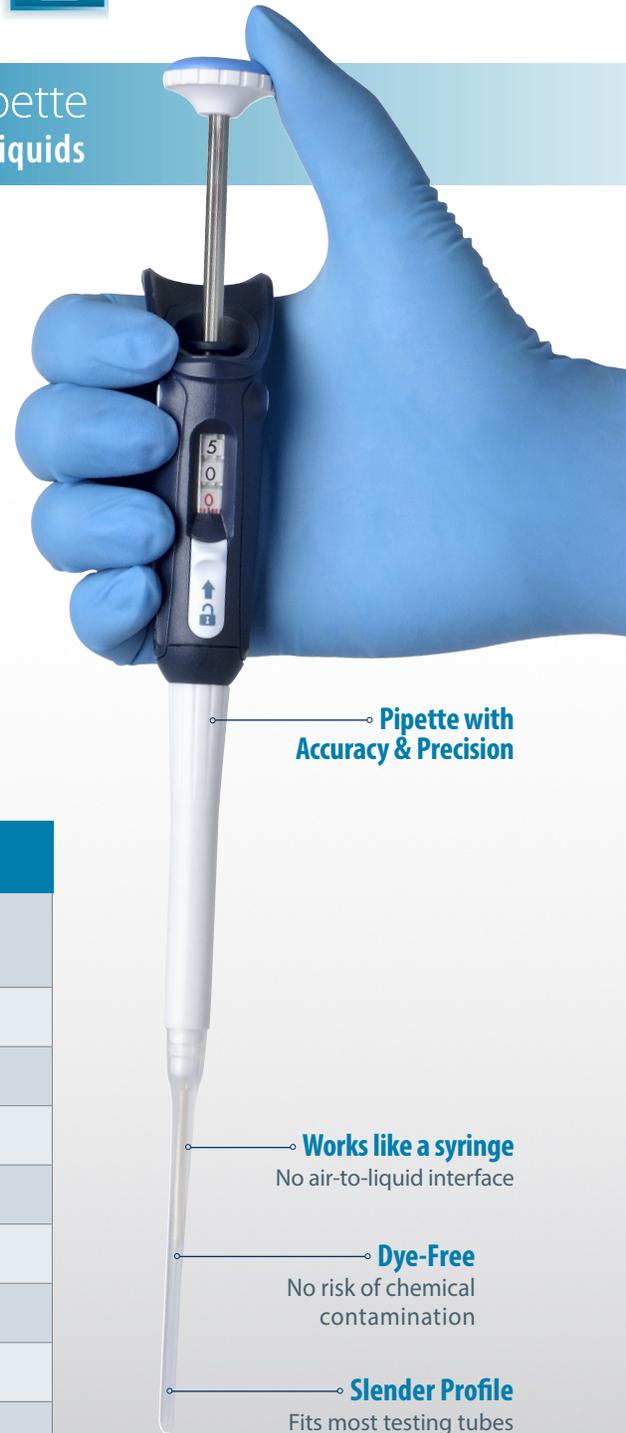

Pipette with Accuracy & Precision

Works like a syringe
No air-to-liquid interface

Dye-Free
No risk of chemical contamination

Slender Profile
Fits most testing tubes

BN1204143



What's New?

New **QuickSnap tip fitting** for easy & instant tip application! MICROMAN® E is the only positive displacement pipette as intuitive and simple to fit as regular pipettes and tips.

New **Repositioned volumeter** — an enhancement from classic MICROMAN®, volume is shown on the front of the pipette for easy view while aspirating and dispensing.

New **Volume-Securing Switch** allows for easy volume adjustment and minimizes the risk of accidentally changing volume during pipetting!

New **Improved Handle Shape** with extended hook allows pipette to rest in the hand comfortably and securely.

New **Dual function push-button** doubles as the tip ejector which prevents user contact with contaminated capillary pistons and allows for easy tip disposal.

Knowing the Difference

Air-displacement pipettes



Three things to remember:

1. Recommended for aqueous samples and for general laboratory work.
2. Always have a cushion of air (dead volume) between the pipette piston and the liquid sample.
3. The piston is a permanent part of the pipette.

Positive-displacement pipettes

Three things to remember:

1. Recommended for problem samples (viscous, dense, volatile, radioactive, corrosive).
2. Direct contact of the piston with the sample (no air cushion).
3. Disposable piston (not permanent part of pipette).

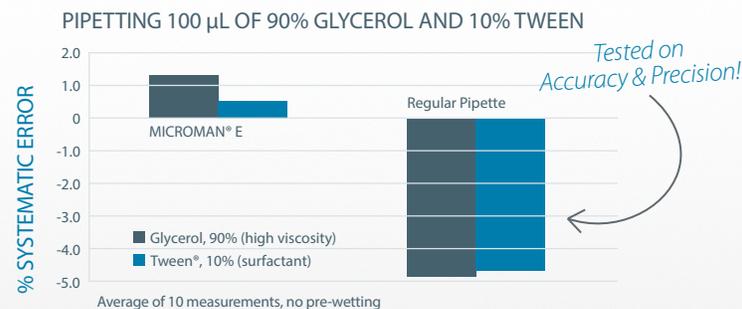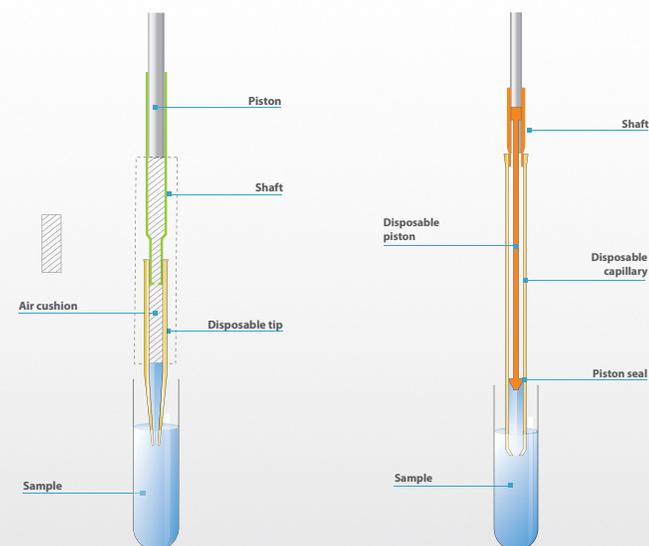

BN1204143

What to Pipette?

With MICROMAN® E pipette everything you can with a standard pipette, PLUS:

- Oils
- Blood
- Glycerol
- Cosmetics
- DNA/RNA
- Radioactive Samples



© Gilson Inc. All rights reserved.

In the United States:

For customer service, call 1-800-766-7000
To fax an order, use 1-800-926-1166
To order online: www.fishersci.com



A Thermo Fisher Scientific Brand